

Getting started with the new MultiLab[®] Pro IDS

1 Connect the power supply




2 Switch on the device




3 Click on "System" & then "Setting"



4 Adapt your language & date/time

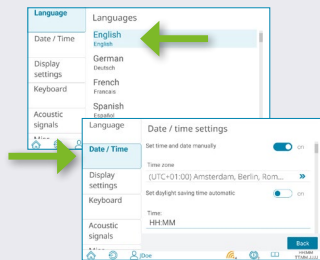

5a Read the manual onboard

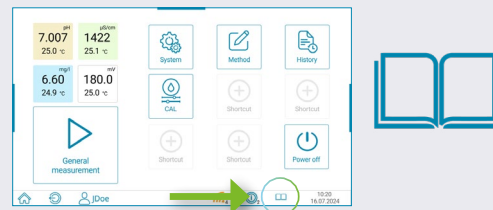


5b Read the manual online




6 Plug-in your sensors



7 Connect your sensors

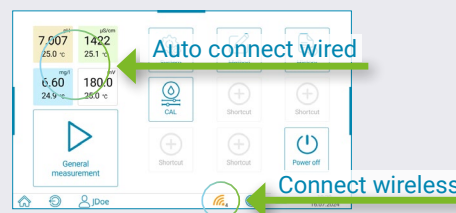

8 Start the measurement

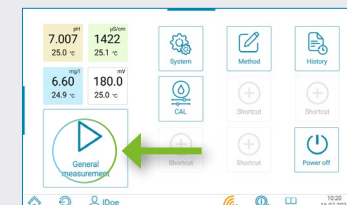

All-in-one solution – measure all electrochemical parameters



pH/ORP

pH & ORP electrodes

- High performance lab or universal electrodes
- Glass or plastic shaft
- Optional built-in temperature sensor
- Penetration / surface / micro / split ring electrodes
- Gel, polymer or liquid filling
- Analog/digital IDS adapter available



BOD

Biochemical Oxygen Demand (BOD) sensors

- Electrochemical or optical versions
- Perfect for BOD₅ testing
- Built-in stirrer
- Compatible with 300 ml BOD bottles



DO

Oxygen sensors


- High performance optical D.O. measurement
- For lab and field applications
- Optical D.O. measurement according to DIN ISO 17289



ISE

Ion-selective electrodes

- Combined ISE & GSE electrodes
- Glass electrodes
- Matrix electrodes
- Solid state electrodes
- Analog/digital IDS adapter available

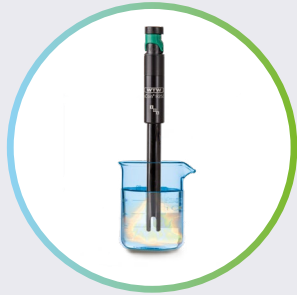


COND

Conductivity cells

- Two-electrode cells for pure and ultrapure water
- Four-electrode cells for general applications
- Graphite or stainless steel electrodes

Check out now our full portfolio of electrodes, sensors and measurement cells:

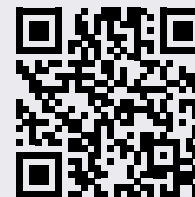

xylem

YSI Incorporated
1725 Brannum Lane
Yellow Springs, OH 45387

Tel +1 937.767.7241
info@ysi.com
YSI.com/xylem-lab-solutions

XA00407 1125

© 2025 Xylem Inc. or its affiliate. All rights reserved.