

Food & Beverage Series

SIMULTANEOUS MEASUREMENT OF GLUCOSE AND SUCROSE IN CORN AND PEAS



Introduction

Glucose (D-glucose) and sucrose concentrations in complex matrices such as corn and peas can be measured directly and quickly using a [YSI 2900 Series Biochemistry Analyzer](#). YSI's unique enzyme technology provides for rapid glucose and sucrose measurements. Measurements are virtually unaffected by color, turbidity, density, pH, or the presence of reducing substances.

When a YSI Biochemistry Analyzer is equipped with a glucose and a sucrose membrane, simultaneous measurement of both analytes is possible. Because glucose interferes with sucrose analysis, it is necessary to follow this protocol when analyzing for sucrose in the presence of glucose.

When a sample is injected into the sample chamber, the sucrose diffuses to the sucrose membrane, which contains invertase, mutarotase, and glucose oxidase. The sucrose is hydrolyzed to α -D-glucose and fructose. The mutarotase allows for the quick equilibrium of glucose between its α and β forms. In the presence of glucose oxidase, the

β -D-glucose (glucose) is oxidized to hydrogen peroxide and D-glucono- δ -lactone. The hydrogen peroxide is detected amperometrically at the platinum electrode surface. The glucose in the sample diffuses to both the glucose and sucrose membranes, which contain glucose oxidase, and is oxidized as well. Subtracting the response of the glucose membrane from the response of the sucrose membrane yields the response due to sucrose alone. The glucose response is taken directly from the glucose membrane. The algorithm in the instrument software calculates the net concentrations. For more information on this system, refer to the Operations Manual.



I. Materials & Setup

- A. YSI Biochemistry Analyzer - equipped with a 2703 Sucrose Membrane, a 2365 Glucose Membrane and 2357 Buffer.
- B. Glucose (2.50 g/L, 25.00 g/L) and Sucrose (5.00 g/L, 25.0 g/L) standard solutions.
- C. Buffer Diluent (40 g/L NaH₂PO₄, 10 g/L Na₂HPO₄ in reagent water).
- D. Connect the YSI Biochemistry Analyzer to a suitable power source.
- E. Perform the instrument and membrane daily checks described in the Operations Manual (Section 5).

- F. Volumetric glassware (Class A recommended).
- G. The following instrument setup is recommended:
Sample Size 10 µL

Probe A Parameters

Chemistry	Glucose
Unit	g/L
Calibrator	2.50 g/L
End Point	30 Sec

Probe B Parameters

Chemistry	Sucrose
Unit	g/L
Calibrator	5.00 g/L
End Point	30 Sec

Autocal Parameters

Temperature	1°C
Time	30 Min
Sample	5 Samples
Cal Shift	2%

II. Method

- A. Weigh up to 25.00 grams of corn or peas to be analyzed. Add about 50 mL buffer diluent.
- B. Transfer the solution to a clean dry blender. Blend until the sample is liquid.
- C. Transfer the sample to a 100 mL volumetric flask using buffer diluent to rinse and dilute. Fill the flask to the mark with buffer and mix. Allow the solution to equilibrate for about twenty minutes before analysis.
- D. Calibrate the YSI Biochemistry Analyzer with a 2.50 g/L glucose and 5.00 g/L sucrose standard solutions.

- E. Check the linearity of the membranes at least once a day by injection of glucose (25.00 g/L) and sucrose (25.0 g/L) linearity check solutions. Refer to the Operations Manual (Section 5) for specifications.
- F. Assay the sample prepared in C by aspiration into the 2900 Series instrument.
- G. Calibrate frequently as described in the Operations Manual (Section 7).

* The linear range of the system is 0 to 25.0 g/L for both glucose and sucrose. The combined concentration of glucose + sucrose cannot exceed 25 g/L. If the sum of the values reported exceeds this, further dilution of the sample is required.

III. Calculations

To calculate % glucose and sucrose, multiply the values reported by the appropriate dilution factor.

Example: 25.10 grams of peas were diluted to 100 mL in a Class A volumetric flask. When assayed, the values reported were 0.128 g/L glucose and 6.91 g/L sucrose.

$$\begin{aligned} \text{\% Glucose:} &= 0.0005 \text{ g glucose / g peas} \\ 1.28 \text{ g/L} \times 0.100 \text{ L} & \\ / 25.10 \text{ g} &= 0.05\% \text{ (w/w)} \end{aligned}$$

$$\begin{aligned} \text{\% Glucose:} &= 0.0275 \text{ g sucrose / g peas} \\ 6.91 \text{ g/L} \times 0.100 \text{ L} & \\ / 25.10 \text{ g} &= 2.75\% \text{ (w/w)} \end{aligned}$$


Ordering Information

YSI Part Numbers:

2900	Series Biochemistry Analyzer
2365	Glucose Membrane Kit
2776	Glucose Standard Solution (2.50 g/L)
2777	Glucose Standard Solution (25.00 g/L)
2703	Sucrose Membrane Kit
2780	Sucrose Standard Solution (5.00 g/L)
2778	Sucrose Standard Solution (25.0 g/L)
2357	Buffer Kit
2363	Potassium Ferrocyanide Test Solution
2392	NaCl Solution (for membrane installation)



APPLICATION NOTE 213LS-02

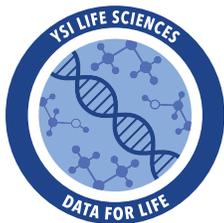

YSI Life Sciences develops and manufactures scientific instruments, sensors and systems that serve a variety of scientific and industrial markets worldwide. YSI has a long history in the life sciences and bioanalytical markets, most notably with our introduction of the world's first commercial whole blood glucose analyzer in 1975. Today there are over 10,000 YSI instruments installed around the world, trusted in critical situations to provide the most accurate data in the shortest time.

YSI, a Xylem brand
1725 Brannum Lane
Yellow Springs, OH 45387

 +1.937.767.7241
 xylem-lab@xylem-inc.com
 YSI.com



[YSI.com/LifeSciences](https://www.ysi.com/LifeSciences)