


# YSI Model 60 Assembly Drawing

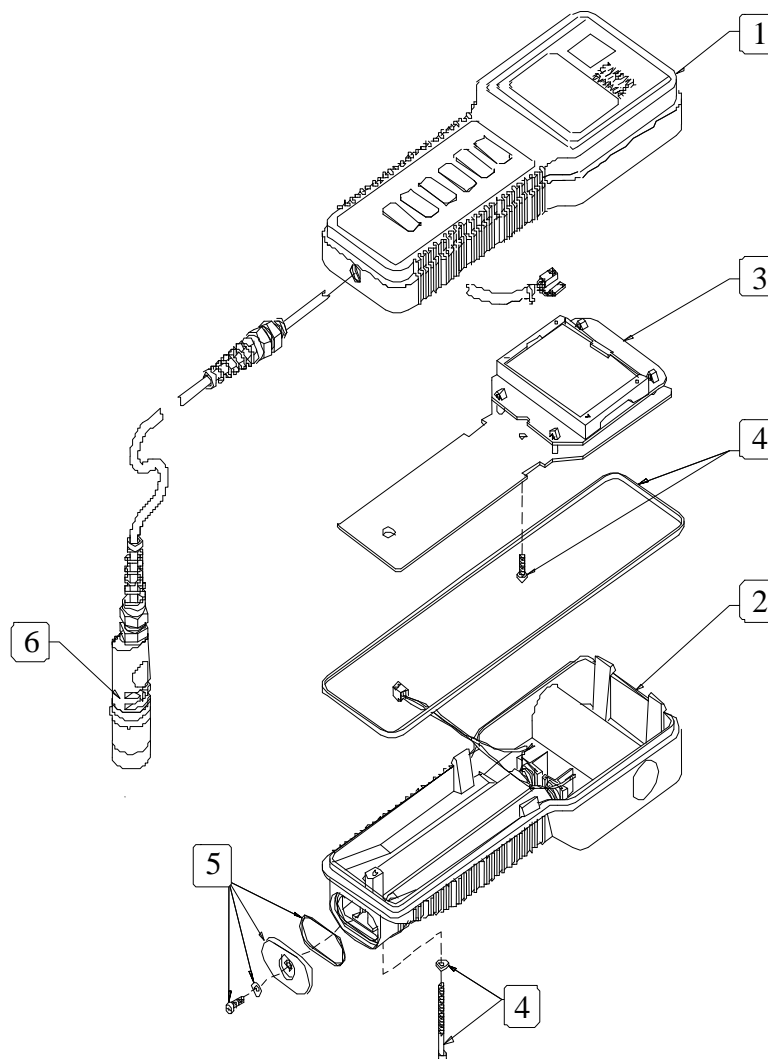

Bubble	YSI Order Number	Description
1	031131	Front Case Assy, W/Keypad
1	055213	Keypad Board Assembly
1	055216	Keypad
2	055242	Rear Case Assy
2	055219	Sponge
3	031128	Board Assy, Main, W/Display
3	055214	LCD Display <u>ONLY</u>
3	115524	LCD Display Kit (includes: LCD, Bracket, Zebra Strips 2 ea.)
4	055204	Kit, Case Hardware
5	055244	Kit, Battery Hardware
6	113135	Cable Assembly, 10 feet
6	113141	Cable Assembly, 25 feet
6	113151	Cable Assembly, 50 feet
6	113171	Cable Assembly, 100 feet
6	131133	Replacement pH Electrode (not shown)
6	031189	100 mL Graduated Cylinder (not shown)