
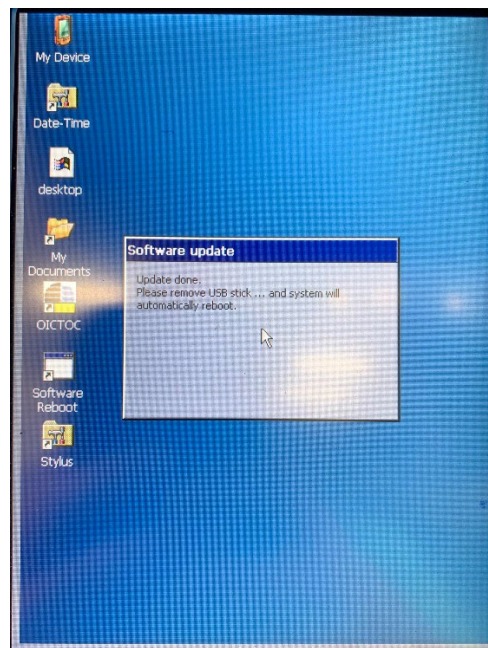


Installation Notes for the TOC 1030 - V1.4.2 FW Patch

Please follow these steps to perform the upgrade:

- **Prepare the USB stick**
 - Download the Zip file for the FW Patch
 - Copy the Patch zip file to a USB stick.
 - Unzip the zip file
 - This will create an **Upgrade** folder on the root folder of the USB stick
- **Prepare the 1030 Instrument**
 - Power OFF the TOC1030 device
 - Remove cover from right side of the 1030 unit
 - Remove the USB stick currently installed in the SOM board
 - Insert the Upgrade USB stick
 - Power ON the TOC1030 device
- **Perform the Upgrade**
 - When the 1030 unit boots up, the desktop screen will appear, and the Upgrade program will launch automatically
 - The upgrade process should take 10-15 seconds.

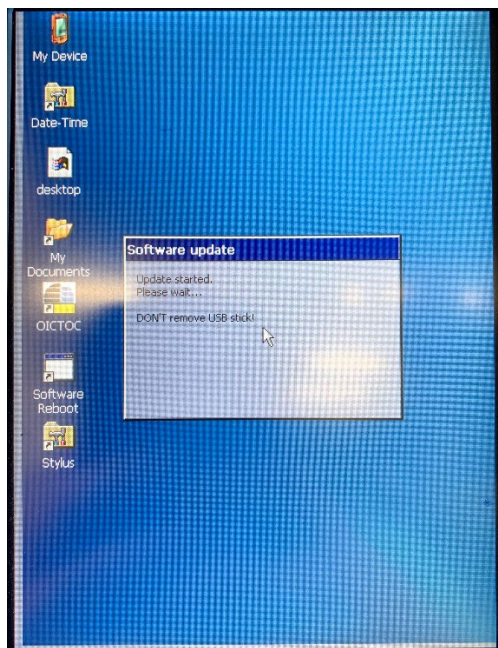

- Once completed, simply remove the USB stick and the 1030 Unit will automatically reboot
- **Reinstall the cover on the 1030 unit**

Update the 1030 Date and Time

- From the **Config-System** menu, click on the **Date/Time** screen
- Edit the Date /Time as needed, then click **Save**